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BEDIENKOMFORT

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USER-FRIENDLINESS

- Rapid production of the cast
- Degassing and cross-linking in a single step
- Simple to use

SAFETY

- High dimensional accuracy
- Bubble-free
- Reproducible and homogenous thickness of the manufactured parts

OPTION

The application of the ZentriForm insert is not limited to silicon wafers. For small batch series or 3D structures customized steel inserts can be used.

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*Microfluidic Chips
Accessories required*

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Centrifuge
ROTOFIX 46, with
heating up to 90 °C
Cat. No. on request

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Add. accessories
Disc spacer in
4 different sizes
Demoulding tool
Cat. No. on request

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Optional accessory
Spin coater
for 4" wafers
Cat. No. 4690

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On request
customized steel
inserts can be
supplied for
3D structures

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www.hettichlab.com
zentriform@hettichlab.com

RAPID PROTOTYPING FOR MICROFLUIDICS

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*The Hettich ZentriForm enables microstructures to be derived **rapidly, simply** and **reliably**. The removal of air bubbles in particular makes the manual production of such structures with silicone a complex process that is susceptible to errors. With the ZentriForm system air bubbles are reliably removed through the centrifugation. The procedure also saves valuable time since curing and degassing are carried out at the same time. The pressure of the silicon against the insert throughout the entire curing process further insures a high dimensional accuracy.*

High reproducibility and a uniform thickness of the parts are the visible result!

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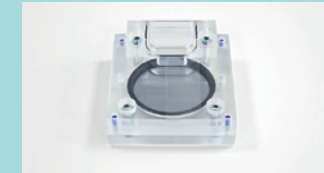
Procedure – rapid, simple and reliable

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


Mount the ZentriForm insert: Insert the master and O-ring and screw the mould together

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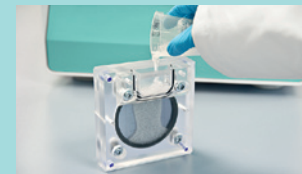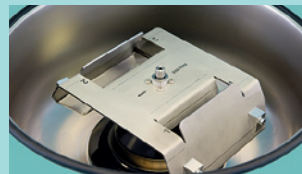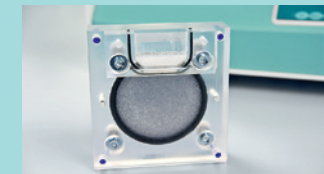Stir the silicon and fill it into the inserts

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Load the rotor and start centrifugation

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Once centrifugation is complete, allow the mould to cool, unscrew it and remove the finished part

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Andreas Hettich GmbH & Co. KG